

How to use GLX For Lab #13

1. Plug a thermostat into a Temperature sensor.
2. Plug four sensors to GLX and turn it on (pressure and temperature will be indicated).
3. Push "Home" button.
4. Select "Sensors" (right bottom).
5. Change "Sample Rate Unit" to "minutes" by pressing check mark and scroll down.
6. Change "Sample Rate" to "5".
7. Since you don't need to see pressure, change it to "Not Visible".
8. Repeat the steps 5 – 7 for all the sensors
9. Press "Home" button to go back to the main menu.
10. Press "Settings" and go to "Auto Power Off".
11. Change it to "60 min".
12. Go back to the main menu and select "Table".
13. Press "F4" and select "Four Columns".
14. Press the play button to start recording.